


Inverted Metallurgical Microscope

GX53

Advanced Microscopy Solutions for Metallurgy Inspection

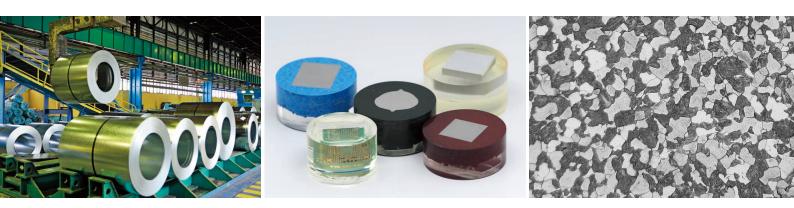





Quick analysis for large or thick sample materials





The GX53 inverted microscope is used for a wide range of applications often seen in the steel, automotive, electronics, and other manufacturing industries. The microscope enables users to inspect polished metals and cross-section samples simply by placing them upside down on the stage. The sample does not need to be leveled and can be thick, large, or heavy.

The GX53 delivers crisp images that can be difficult to capture using conventional microscopy observation methods. When combined with OLYMPUS Stream image analysis software, the microscope streamlines the inspection process from observation to image analysis and reporting.

Streamline Your Inspection Process

Fast Inspections, Advanced Functionality

Quickly observe, measure, and analyze metallurgical structures.

User-Friendly

Even novice operators can comfortably make observations, analyze results, and create reports.

Advanced Imaging Technology

Our proven optics and imaging technology deliver clear images and reliable results.

Modular

Choose the components you need for your application.


Fast Inspections, Advanced Functionality

Advanced Analysis Tools

The GX53 microscope's various observation capabilities provide clear, sharp images, so users can reliably detect defects in their samples. OLYMPUS Stream image analysis software's new illumination techniques and image acquisition options give users more choices for evaluating their samples and documenting their findings.

The Invisible Becomes Visible: MIX Technology


MIX technology produces unique observation images by combining darkfield with another observation method, such as brightfield or polarization. MIX observation enables users to view samples that are difficult to see with conventional microscopes, and represents even small height differences of sample surfaces. The circular LED illuminator used for darkfield observation has a directional darkfield function where one or more quadrants are illuminated at a given time. This reduces a sample's halation and is useful for visualizing its surface texture.

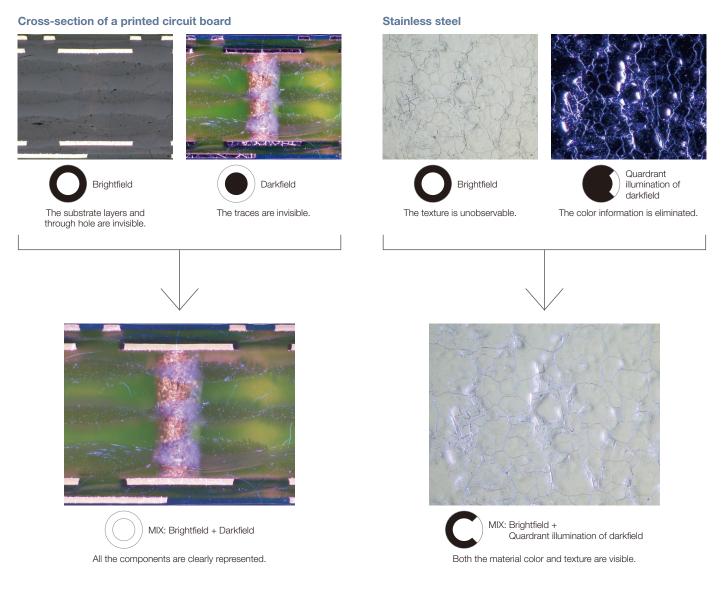

Easily Create Panoramic Images: ST Instant MIA

With multiple image alignment (MIA), users can stitch images together quickly and simply by moving the XY knobs on the manual stage-a motorized stage is optional. OLYMPUS Stream software uses pattern recognition to generate a panoramic image, which is suitable for inspections of carburizing and metal-flow conditions.

Metal flow of a bolt

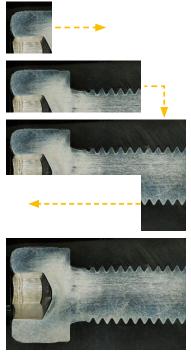

The full condition of metal flow can be seen.

Create All-in-Focus Images: EFI

OLYMPUS Stream software's extended focus imaging (EFI) function captures images of samples whose height extends beyond the depth of focus. EFI stacks these images together to create a single all-in-focus image of the sample. Even when analyzing a cross-section sample with an uneven surface, EFI creates fully-focused images. EFI works with either a manual or motorized Z-axis and creates a height map to easily visualize structures.

Resin parts






Fully focused image

Capture Both Bright and Dark Areas Using HDR

Using advanced image processing, high dynamic range (HDR) adjusts for differences in brightness within an image to reduce glare. It also helps boost the contrast in low-contrast images. HDR can be used to observe minute structures in electric devices and identify metallic grain boundaries.

Gold plate

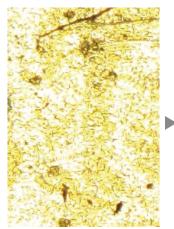

Some areas have glare.



Both dark and bright areas are clearly exposed using HDR.

Chromium diffusion coating



Low contrast and unclear.




Enhanced contrast with HDR.

OLYMPUS Stream Software – Optimized for Materials Science

Material inspection, measurement, and analysis are required to comply with industrial standards as well as internal operating procedures. Together, the GX53 microscope and OLYMPUS Stream software support metallurgical analysis methods that comply with different industrial standards. With step-by-step operator guidance, users can analyze their samples quickly and easily.

Particle Analysis – Count and Measure Solution

Detecting objects and measuring size distribution are among the most important applications in digital imaging. OLYMPUS Stream software's Count and Measure solution uses advanced threshold methods to reliably separate objects, such as particles and scratches, from the background. More than 50 different object measurement and classification parameters are available including shape, size, position, and pixel properties.

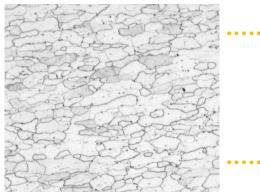


Etched steel microstructure (original image)

Conventional software

Unclear grain boundaries

OLYMPUS Stream

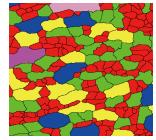

Grain boundaries are clearly detected




Grain classification results


Grain Sizing in a Microstructure

Users can measure the grain size and analyze the microstructure of aluminum, steel crystal structures, such as ferrite and austenite, and other metals. Supported standards: ISO, GOST, ASTM, DIN, JIS, GB/T

Microstructure of ferritic grains

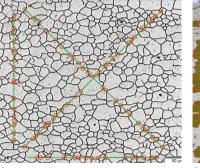

Grain sizing intercept solution

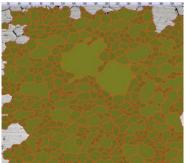

Grain sizing planimetric solution

Evaluating Graphite Nodularity

The software can be used to evaluate the graphite nodularity and content in cast iron samples (nodular and vermicular). The form, distribution, and size of graphite nodes can be classified.

Supported standards: ISO, NF, ASTM, KS, JIS, GB/T

Ductile cast iron showing nodular graphite




Cast iron solution

Rating Nonmetallic Inclusion Content in High-Purity Steel

Classify nonmetallic inclusions using an captured image of the worst field or inclusion that you have manually located on the sample.

Supported standards: ISO, EN, ASTM, DIN, JIS, GB/T, UNI

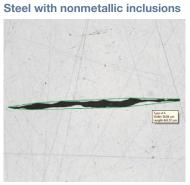


Inclusion worst field solution

Compare Images of Your Sample and Reference Images

Easily compare live or still images with auto-scaled reference images. This solution includes reference images in accordance with various standards. The solution also supports multiple modes, including live overlay display and side-by-side comparison. Additional reference images can be purchased separately. Supported standards: ISO, EN, ASTM, DIN, SEP

Steel with nonmetallic inclusions

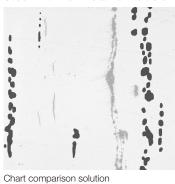

Microstructure with ferritic grains



Chart comparison solution

Material Solution Specifications*

| Solutions | Supported standards |
|-----------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Grain intercept | ISO 643: 2012, JIS G 0551: 2013, JIS G 0552: 1998, ASTM E112: 2013, DIN 50601: 1985, GOST 5639: 1982, GB/T 6394: 2002 |
| Grain planimetric | ISO 643: 2012, JIS G 0551: 2013, JIS G 0552: 1998, ASTM E112: 2013, DIN 50601: 1985, GOST 5639: 1982, GB/T 6394: 2002 |
| Cast iron | ISO 945-1: 2010, ISO 16112: 2017, JIS G 5502: 2001, JIS G 5505: 2013, ASTM A247: 16a, ASTM E2567: 16a, NF A04-197: 2004, GB/T 9441: 2009, KS D 4302: 2006 |
| Inclusion worst field | ISO 4967 (method A): 2013, JIS G 0555 (method A): 2003, ASTM E45 (method A): 2013, EN 10247 (methods P and M): 2007, DIN 50602 (method M): 1985, GB/T 10561 (method A): 2005, UNI 3244 (method M): 1980 |
| Chart comparison | ISO 643: 1983, ISO 643: 2012, ISO 945: 2008, ASTM E 112: 2004, EN 10247: 2007, DIN 50602: 1985, ISO 4505: 1978, SEP 1572: 1971, SEP 1520: 1998 |
| Coating thickness | EN 1071: 2002, VDI 3824: 2001 |

* Please see the OLYMPUS Stream brochure for more detailed information.

User-Friendly

A Design That Emphasizes User Comfort

The microscope's ergonomic design helps users stay comfortable while they work, contributing to a more efficient inspection. When used with OLYMPUS Stream software, operators can easily acquire images of diverse samples, conduct a variety of analyses, and generate professional reports.

Maintain a comfortable posture

The tilting observation tube's extensive range and adjustable eyepoint enable operators to sit or stand at the microscope in a comfortable posture.



Observe large, heavy samples

Samples up to 5 kg can be inspected simply by placing the polished surface on the stage.

Helps prevent objective collisions

The stage mirror helps make it easier to adjust the observation point and objective magnification. It also helps prevent the objective from colliding with the sample.



Easily switch observation methods

The microscope supports brightfield, darkfield, differential interference contrast (DIC), and simple polarized light observations. Use a dedicated level to quickly switch between brightfield and darkfield.



Add DIC simply by adding a slider.

Instantly record observation images

With the touch of a button (optional), observed images can be instantly saved.



Convenient hand switch

Control MIX illumination, the objectives, and OLYMPUS Stream functions using the available hand switch.




Easily control the stage during observation

Use the dedicated handle to control the stage while you are looking through the eyepieces.




Easily Restore Microscope Settings: Coded Hardware

Coded functions integrate the microscope's hardware settings with OLYMPUS Stream image analysis software. The observation method, illumination intensity, and magnification can be recorded by the software and stored with the associated images. Since the settings can easily be reproduced, different operators can conduct the same quality inspections with limited training.

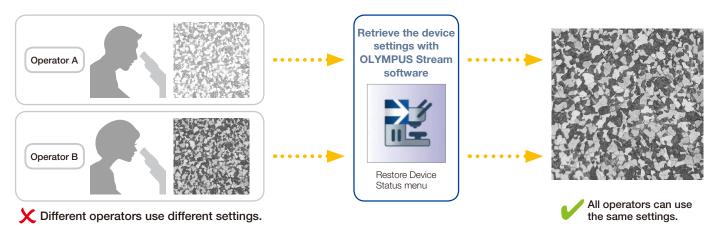

User Guidance Helps Simplify Advanced Analysis


The software guides users step-by-step through an inspection process that complies with the chosen industrial standard. Operators at any experience level can quickly and easily conduct advanced analysis by following the on-screen guidance.

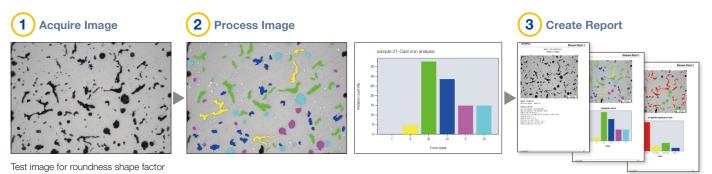

Efficient Report Generation

Creating a report can often take longer than capturing the image and taking the measurements. OLYMPUS Stream software provides intuitive report creation to repeatedly produce smart and sophisticated reports based on predefined templates. The software can be configured so that the magnification is printed along with individual images.



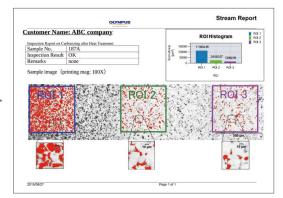

8


Advanced Imaging Technology


Proven optics and digital imaging technology deliver quality inspection data

Olympus' history of developing high-quality optics and advanced imaging capabilities has led to quality microscopes that offer exceptional measurement accuracy.

Reliable Optical Performance: Wavefront Aberration Control


The optical performance of objective lenses directly impacts the quality of the observation images and analysis results. Olympus UIS2 high-magnification objectives are designed to minimize wavefront aberrations, delivering reliable optical performance.

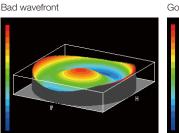

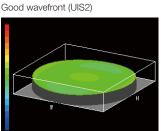

Consistent Color Temperature: High-Intensity White LED Illumination

The GX53 microscope utilizes a high-intensity white LED light source for reflected and transmitted illumination. The LED maintains a consistent color temperature regardless of intensity for reliable image quality and color reproduction. The LED system provides efficient, long-life illumination that is ideal for materials science applications.

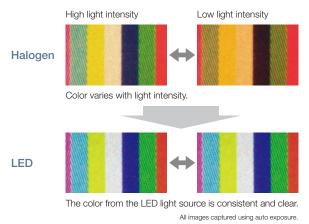


Entirely Clear Image: Image Shading Correction




OLYMPUS Stream software features shading correction to mitigate shading in the corners of an image. When used with intensity threshold settings, shading correction provides a more precise analysis.

Stainless steel (Binarized image)

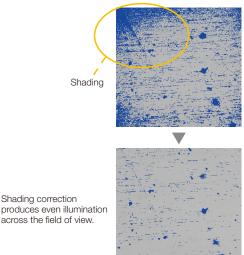

Precise Measurements: Auto Calibration

Similar to digital microscopes, automatic calibration is available when using OLYMPUS Stream software. Auto calibration helps eliminate the impact of human variability on the calibration process, leading to more reliable measurements. The software automatically calculates the correct calibration from an average of multiple measurement points, minimizing variance and

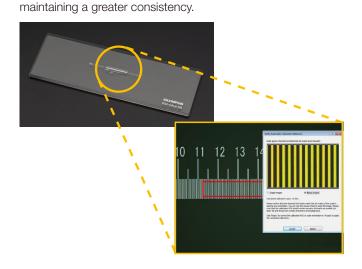


Applications

Reflected light microscopy spans a range of applications and industries. Below are just a few examples of what can be achieved using different observation methods.

Polished sample of AlSi



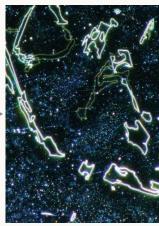


Brightfield

Darkfield

Brightfield is a common observation method to observe reflected light from a sample by illuminating it straight on.

Darkfield is used to observe scattered or diffracted light from a sample, so imperfections clearly stand out. Inspectors can identify even minute scratches or flaws.

Spheroidal graphite cast iron

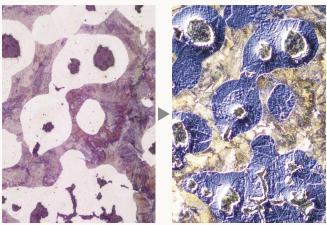

Brightfield


DIC observation

Differential interference contrast (DIC) is an observation technique where the height of a sample, normally not detectable in brightfield, is visible as a relief, similar to a 3D image with improved contrast. It is ideal for inspections of samples that have very minute height differences, including metallurgical structures and minerals.

Aluminum alloy




Brightfield

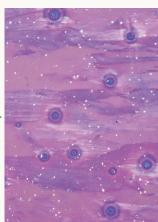

Polarized light observation

Polarized light observation represents a material's texture and crystal condition brightly. It is suitable for metallurgical structures such as the growth pattern of graphite on nodular cast iron and minerals.

Electronic device

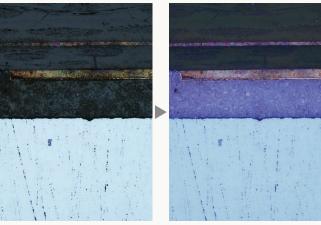

Brightfield

MIX observation: Brightfield + Darkfield

MIX observation combines brightfield and darkfield illumination methods, showing both the sample's color and structure.

The above MIX observation image clearly reproduces the device's color and texture as well as the condition of the adhesive layer.

Customizable

Choose the Components You Need

The GX53 microscope is designed to enable users to choose a variety of optical components to suit individual inspection and application requirements. The system can utilize all available observation methods. Users can also select from a variety of OLYMPUS Stream image analysis packages to meet image acquisition and analysis needs.

GX53 Reflected/Transmitted Light Combination

The GX53 microscope frame can be configured for both reflected and transmitted light with manual, coded, or motorized components.



Scales for Metallurgical Analysis

Glass scales can be inserted into the eyepiece to conduct observations that comply with industry standards. Grain size reticles, squared circles, and calibration scales are also available for each objective.

Scale slider

| 1 | GX-SLM | Scale slider, attachable 3 glass scales maximum |
|---|--------------|-------------------------------------------------------------------------------------------------------------|
| 2 | GX51-SLMG5 | Scale glass for 5× objective, scale length: 200 µm |
| 3 | GX51-SLMG10 | Scale glass for 10× objective, scale length: 100 µm |
| 4 | GX51-SLMG20 | Scale glass for 20× objective, scale length: 50 µm |
| 5 | GX51-SLMG50 | Scale glass for 50× objective, scale length: 10 µm |
| 6 | GX51-SLMG100 | Scale glass for 100×, scale length: 10 µm |
| 7 | GX51-SLMGS | Grain size scale, applied to JIS G 0551, ISO 643 and ASTM E112 AUSTENITE GRAINS IN STEEL PLATE IV No.1 to 8 |
| 8 | GX51-SLMGH | Lattice pattern, applied to JIS G 0555 |
| 9 | GX-SLMG | Parfocal glass to adjust the light path length |





Build Your System Your Way

Microscope Frame

The GX53 microscope has a built-in power supply for reflected light. The camera adaptor port at the front of the microscope enables users to display live and captured images without using a trinocular tube. Choose various accessories such as a stage mirror that enables users to check the observation position and the magnification of the objectives.

Microscope frames

| : Possible | | Reflected light | Transmitted light | | | | | | |
|------------|-------------|-----------------------------------|-------------------|--|--|--|--|--|--|
| 1 | GX53F | | | | | | | | |
| Access | Accessories | | | | | | | | |
| 2 | CK40M-MS | Observation position check mirror | | | | | | | |
| - | COVFR-021 | Dust cover for GX53 system | | | | | | | |

Transmitted Illumination Unit

Condensers collect and focus transmitted light and are used for transmitted light observation.

| 1 | IX2-ILL100 | Stand for transmitted illumination, attachable BF/POL lamp housing fo LED (BX3M-LEDT) and halogen (U-LH100L-3) | | | | |
|---|------------|-------------------------------------------------------------------------------------------------------------------|--|--|--|--|
| 2 | PMG3-LWCD | Condenser for transmitted light observation, condenser (NA 0.6, WD 12 mm) with aperture stop | | | | |

Light Sources

Choose the light source and power supply you need to illuminate your sample. Choose the appropriate light source for your observation method.

Standard LED light source configuration

| 1 | BX3M-LEDR | LED lamp housing for reflected light |
|---|------------|-------------------------------------------------------------------------|
| 2 | BX3M-LEDT | LED lamp housing for transmitted light |
| 3 | BX3M-PSLED | Power supply for LED lamp housing (required for transmitted light only) |

High intensity light source configuration

| | - | | | | | |
|-------|-----------|----------------|----------------------------------------------------------------------------------------|--|--|--|
| 4 | MX-HGAD | | High intensity light adapter | | | |
| 5 | LU-I | LGAD | Liquid light guide adapter | | | |
| 6, 7 | L | J-LLG150 (300) | Liquid light guide, length: 1.5 m (3 m) | | | |
| 8 | -U-HGLGPS | | High intensity light source (mercury), including one SHI-130OL in the standard package | | | |
| - | | LSHI-130OL | 130 W mercury lamp | | | |
| 9, 10 | LU-L | H100HG (HGAPO) | Mercury lamp housing, chromatic aberration correction type | | | |
| - | | -USH-103OL | 100 W mercury lamp | | | |
| 11 | LU-CLA | | Flexible extension handle for mercury lamp housing | | | |
| 12 | -U-RFL-T | | Power supply for 100 W mercury lamp | | | |
| 13 | L | J-CST | Optical axis adjustment sample for mercury lamp housing | | | |

Halogen light source configuration

| 14 | U-LH100L-3 | Halogen lamp housing | | | | |
|--------|---------------------|-------------------------------------------------------------------------------------------------------|--|--|--|--|
| - | -12V100W HAL (-L) | 100 W halogen lamp (long life type) | | | | |
| 15 | L _{U-RMT} | Extender cable for halogen lamp housing, cable length 1.7 m (requires cable extension when necessary) | | | | |
| 16, 17 | | 100 V (200 V) specification power supply for 100 W/50 W halogen lamp | | | | |
| 18 | L _{TH4-HS} | Hand switch to change the light intensity of halogen (dimmer TH4-100 (200) without hand switch) | | | | |

Double lamp housing configuration

| 19 | Ų-DULHA | Dual lamp housing attachment | | | | |
|----|-----------------------------------------------------------------------------------------------------|------------------------------|--|--|--|--|
| | High intensity light source configuration (MX-HGAD is not required when using U-LH100HG (HGAPO)) | | | | | |
| | BX3M-LEDR (with standard LED light source configuration) | | | | | |
| | Halogen light source configuration | | | | | |







Tubes

Select tubes for imaging through the eyepieces or for use with a camera. Choose the tube you need by imaging type and level of ergonomic comfort.

| | | FN (mm) | Туре | Angle type | Image | Diopter adjustment mechanism | Turret mechanism |
|---|-----------|---------|------------|------------|---------|------------------------------------|---------------------|
| 1 | U-BI90 | 22 | Binocular | Fixing | Reverse | Right only | - |
| 2 | U-BI90CT | 22 | Binocular | Fixing | Reverse | Right only | 4 positon* |
| 3 | U-TBI90 | 22 | Binocular | Tilting | Reverse | Right only | _ |
| 4 | U-TR30H-2 | 22 | Trinocular | Fixing | Reverse | Right only | - |

*4 positions are O, CT, O, and S.

(O: Empty, CT: Centering telescope for adjustment of aperture stop, S: Shutter preventing light from eyepiece.)

Eyepieces

Eyepiece for viewing directly into the microscope. Select based on the desired field of view.

| : Possible | | FN (mm) | Diopter adjustment mechanism | Built-in cross reticle |
|------------|--------------|------------|---------------------------------|------------------------|
| 1 | WHN10X | 22 | | |
| 2 | WHN10X-H | 22 | | |
| 3 | CROSS WHN10X | 22 | | |

Intermediate Tubes

Various accessories for multiple purposes. For use between the tube and microscope frame.

| 1 | U-CA | Magnification changer (1×, 1.25×, 1.6×, 2×) | | | | |
|---|--------|---------------------------------------------|--|--|--|--|
| 2 | U-ECA | Agnification changer (1×, 2×) | | | | |
| 3 | U-EPA2 | Eyepoint adjuster : + 30 mm | | | | |
| 4 | GX-SPU | Attachable camera adapter with side port | | | | |
| 5 | IX-ATU | Attachable tube : U-TR30H-2 | | | | |

Camera Adapters


Adaptors are used to add a camera. Select the adaptor based on the field of view and magnification. The actual observation range can be calculated using the following formula: actual field of view (diagonal mm) – viewing field (viewing number) / objective magnification.

| | | Magnification | Centering adjustment | adjustment (field | | mera image area ld number) (mm) | | |
|-----------|-----------------------------|---------------|----------------------|-------------------|-----------|------------------------------------|-----------|--|
| | | | (mm) | 2/3 in. | 1/1.8 in. | 1/2 in. | unit | |
| 1 | GX-TV0.7XC | 0.7 | - | 15.3 | 12.6 | 11.4 | GX53F | |
| 2 | GX-TV0.5XC | 0.5 | - | 21.4 | 17.6 | 16 | GX53F | |
| 3 | U-TV1X-2 with U-CMAD3 | 1 | - | 10.7 | 8.8 | 8 | GX-SPU | |
| 4 | U-TV1XC | 1 | ø2 | 10.7 | 8.8 | 8 | GX-SPU | |
| 5 | U-TV0.63XC | 0.63 | - | 17 | 14 | 12.7 | GX-SPU | |
| 6 | U-TV0.5XC-3 | 0.5 | - | 21.4 | 17.6 | 16 | GX-SPU | |
| 7 | U-TV0.35XC-2 | 0.35 | - | - | _ | 22 | GX-SPU | |
| 8 | U-TV0.25XC* | 0.25 | - | - | - | - | GX-SPU | |
| 9, 10, 11 | IX-TVAD with U-FMT/U-CMT | 1 | - | 10.7 | 8.8 | 8 | U-TR30H-2 | |

For information on digital cameras, please visit our website at http://www.olympus-ims.com/en/microscope/dc/ * A camera can be attached when the image area (field number) is less than 1/3 inch.





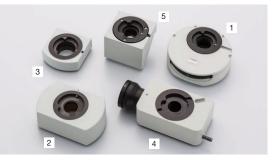



Nosepieces

Nosepieces are used to attach objectives and sliders. Choose your nosepiece based on the number of objectives you want to attach, objective type, and whether or not you are using a slider attachment.

| | : Possible | Туре | Holes | BF | DF | DIC | MIX | ESD | Number of centering holes |
|----|---------------|--------|-------|----|----|-----|-----|-----|---------------------------|
| 1 | U-5RE-2 | Manual | 5 | | | | | | |
| 2 | U-5RES-ESD | Coded | 5 | | | | | | |
| 3 | U-P4RE | Manual | 4 | | | | | | 4 |
| 4 | U-D6RE | Manual | 6 | | | | | | |
| 5 | U-D6RE-ESD-2 | Manual | 6 | | | | | | |
| 6 | U-P6RE | Manual | 6 | | | | | | 2 |
| 7 | U-D7RE | Manual | 7 | | | | | | |
| 8 | U-D6RES | Coded | 6 | | | | | | |
| 9 | U-D7RES | Coded | 7 | | | | | | |
| 10 | U-5BDRE | Manual | 5 | | | | | | |
| 11 | U-D5BDRE | Manual | 5 | | | | | | |
| 12 | U-P5BDRE | Manual | 5 | | | | | | 2 |
| 13 | U-D6BDRE | Manual | 6 | | | | | | |
| 14 | U-D5BDRES-ESD | Coded | 5 | | | | | | |
| 15 | U-D6BDRES-S | Coded | 6 | | | | | | |



Sliders

Select the slider to complement traditional brightfield observation. The DIC slider provides topographic information about the sample with options to maximize contrast or resolution. The MIX slider provides illumination flexibility with a segmented LED source in the darkfield path.

| | Туре | | Amount of shear | Recommended objectives |
|---|----------|------------|-----------------|-------------------------------------------|
| 1 | U-DICR | Standard | Medium | MPLFLN, MPLAPON, LMPLFLN, and LCPLFLN-LCD |
| 2 | U-DICRH | Resolution | Small | MPLFLN, MPLAPON |
| 3 | U-DICRHC | Contrast | Large | LMPLFLN and LCPLFLN-LCD |

MIX slider for MIX observation

| | | Туре | Available objectives |
|---|--------|------------|--------------------------------|
| 4 | U-MIXR | MIX slider | MPLFLN-BD, LMPLFLN-BD, MPLN-BD |

Control Box Hand Switches

Control boxes for connecting the microscope's hardware with a PC and hand switches for hardware display and control.

Control box

| 1 | BX3M-CBFM | Control box for the BXFM system | |
|--------|-----------|-----------------------------------------------------------------------------------------------------------------------------------------------|--|
| 2 | GX-IFRES | Box for OB indicator of the hand switch BX3M-HS; If the GX-IFRES connects to BX3M-CBFM, U-CBS is not needed when using OLYMPUS Stream/DP2-SAL | |
| 3 | U-CBS | Control box for coded functions | |
| Hand s | witch | | |
| 4 | BX3M-HS | MIX observation control, indicator of coded/motorized hardware, programmable software function button of OLYMPUS Stream | |
| 5 | U-HSEXP | Operate a camera's shutter | |
| Cable | | | |
| - | U-MIXRCBL | U-MIXR cable, cable length: 0.5 m | |





Stages

Stages and stage plates for sample placement. Select based on sample shape and size.

Stages

| 1 | IX2-SFR | Flexible right handle stage, the handle grip is about 260 mm below the stage surface | |
|---|----------|----------------------------------------------------------------------------------------------------------------------------------------------|--|
| 2 | GX-SFR | Flexible right handle stage, the handle grip is about 280 mm below the stage surface | |
| 3 | GX-SVR | Right handle stage | |
| 4 | IX2-GS | Gliding stage, the stage plate is incorporated (diameter: ø110 mm, shape of hole: ø25 mm teardrop, material: aluminum alloy) | |
| 5 | IX-SVL-2 | Flexible left handle (short) stage, stage plate is incorporated (diameter: 110 mm, shape of hole: ø25 mm teardrop, material: aluminum alloy) | |

Stage plates

| | | Plate area | Hole type | Material |
|---|------------|------------|------------------------|-------------|
| 6 | CK40-CPG30 | ø110 mm | Diameter ø30 mm | Glass |
| 7 | IX-CP50 | ø110 mm | Diameter ø50 mm | Brass |
| 8 | IX2-GCP | ø110 mm | Teardrop ø25 mm | Brass |
| | GX-CP | a110 mm | Teardrop ø12 mm | Brass |
| 9 | GX-CP | ø110 mm | Long hole (74 × 25 mm) | Amber alloy |



Optical Filters

Optical filters convert sample exposure light to various types of illumination. Select the appropriate filter for your observation requirements.

BF, DF, FL

| 1, 2, 3 | U-25ND50, 25, 6 | Transmittance 50%/25%/6% |
|---------|--------------------|---------------------------------------------------------------|
| 4 | U-25LBD | Daylight color filter |
| 5 | U-25LBA | Halogen color filter |
| 6 | U-25IF550 | Green filter |
| 7 | U-25L42 | UV cut filter |
| 8 | U-25Y48 | Yellow filter |
| 9 | U-25FR | Frost filter |
| 10 | GX-FSL | Used by combining GX51 filters, attachable filter quantity: 3 |
| 11, 12 | -25ND25, 6 | ø25 mm transmittance 25%/6% |
| 13 | -25LBD | ø25 mm daylight color filter |
| 14 | -25IF550 | ø25 mm green filter |
| 15 | L _{25Y48} | ø25 mm yellow filter |

POL, DIC

| , | - | | | | |
|----|---------------------------------------------------------------------|------------------------------------------------------------------------------|--|--|--|
| 16 | GX-AN Analyzer for reflected light; polarization direction is fixed | | | | |
| 17 | GX-AN360 | Analyzer for reflected light; polarization direction is 360 degree rotatable | | | |
| 18 | GX-PO3 | Polarizer for reflected light; polarization direction is fixed | | | |
| 19 | GX-POTP | Tint plate polarizer for reflected light; polarization direction is fixed | | | |
| | | | | | |

Transmitted light

| 20 | U-POT | ø45 mm polarizer filter |
|--------|-------------|----------------------------------------------------|
| 21 | 43IF550-W45 | ø45 mm green filter for transmitted light |
| 22 | 45-LBD-IF | ø45 mm daylight color filter for transmitted light |
| 23, 24 | 45-ND25, 6 | ø45 mm transmittance 25%/6% for transmitted light |
| | · | |

Other

| | 25 | U-25 | Empty filter, used by combining user's ø25 mm filters |
|---|----|------|-------------------------------------------------------|
| _ | | | |



UIS2 Objectives

Objectives magnify the sample. Select the objective that matches the working distance, resolving power, and observation method for the application.

| Objectives | | Magnifications | NA | W.D. (mm) | Cover grass thickness*2 (mm) | Resolution*3 (µm) |
|---------------|----|----------------|------|--------------|------------------------------------|----------------------|
| | 1 | 50X | 0.95 | 0.35 | 0 | 0.35 |
| MPLAPON | 2 | 100X | 0.95 | 0.35 | 0 | 0.35 |
| | 3 | 1.25X*4*5 | 0.04 | 3.5 | 0-0.17 | 8.39 |
| | 4 | 2.5X*5 | 0.08 | 10.7 | 0-0.17 | 4.19 |
| | 5 | 5X | 0.15 | 20 | 0-0.17 | 2.24 |
| | 6 | 10X | 0.30 | 11 | 0-0.17 | 1.12 |
| MPLFLN | 7 | 20X | 0.45 | 3.1 | 0 | 0.75 |
| | 8 | 40X*1 | 0.75 | 0.63 | 0 | 0.45 |
| | 9 | 50X | 0.80 | 1 | 0 | 0.42 |
| | 10 | 100X | 0.90 | 1 | 0 | 0.37 |
| | 11 | 20X | 0.25 | 25 | 0-0.17 | 1.34 |
| SLMPLN | 12 | 50X | 0.35 | 18 | 0 | 0.96 |
| | 13 | 100X | 0.60 | 7.6 | 0 | 0.56 |
| | 14 | 5X | 0.13 | 22.5 | 0-0.17 | 2.58 |
| | 15 | 10X | 0.25 | 21 | 0-0.17 | 1.34 |
| LMPLFLN | 16 | 20X | 0.40 | 12 | 0 | 0.84 |
| | 17 | 50X | 0.50 | 10.6 | 0 | 0.67 |
| | 18 | 100X | 0.80 | 3.4 | 0 | 0.42 |
| | 19 | 5X | 0.10 | 20 | 0-0.17 | 3.36 |
| | 20 | 10X | 0.25 | 10.6 | 0-0.17 | 1.34 |
| MPLN*4 | 21 | 20X | 0.40 | 1.3 | 0 | 0.84 |
| | 22 | 50X | 0.75 | 0.38 | 0 | 0.45 |
| | 23 | 100X | 0.90 | 0.21 | 0 | 0.37 |
| | 24 | 20X | 0.45 | 8.3-7.4 | 0-1.2 | 0.75 |
| LCPLFLN-LCD | 25 | 50X | 0.70 | 3.0-2.2 | 0-1.2 | 0.48 |
| | 26 | 100X | 0.85 | 1.2-0.9 | 0-0.7 | 0.39 |
| | 27 | 5X | 0.15 | 12 | 0-0.17 | 2.24 |
| | 28 | 10X | 0.30 | 6.5 | 0-0.17 | 1.12 |
| | 29 | 20X | 0.45 | 3 | 0 | 0.75 |
| MPLFLN-BD*6 | 30 | 50X | 0.80 | 1 | 0 | 0.42 |
| | 31 | 100X | 0.90 | 1 | 0 | 0.37 |
| | 32 | 150X | 0.90 | 1 | 0 | 0.37 |
| | 33 | 5X | 0.15 | 12 | 0-0.17 | 2.24 |
| | 34 | 10X | 0.25 | 6.5 | 0-0.17 | 1.34 |
| MPLFLN-BDP*6 | 35 | 20X | 0.40 | 3 | 0 | 0.84 |
| | 36 | 50X | 0.75 | 1 | 0 | 0.45 |
| | 37 | 100X | 0.90 | 1 | 0 | 0.37 |
| | 38 | 5X | 0.13 | 15 | 0-0.17 | 2.58 |
| | 39 | 10X | 0.25 | 10 | 0-0.17 | 1.34 |
| LMPLFLN-BD*6 | 40 | 20X | 0.40 | 12 | 0 | 0.84 |
| | 41 | 50X | 0.50 | 10.6 | 0 | 0.67 |
| | 42 | 100X | 0.80 | 3.3 | 0 | 0.42 |
| | 43 | 5X | 0.10 | 12 | 0-0.17 | 3.36 |
| | 44 | 10X | 0.25 | 6.5 | 0-0.17 | 1.34 |
| MPLN-BD*4*6*7 | 45 | 20X | 0.40 | 1.3 | 0 | 0.84 |
| | 46 | 50X | 0.75 | 0.38 | 0 | 0.45 |
| | 47 | 100X | 0.90 | 0.21 | 0 | 0.37 |





 $\star 1$ The MPLFLN40X objective is not compatible with the differential interference contrast microscopy.


*2 0: For viewing specimens without a cover glass.

*3 Resolutions calculated with aperture iris diaphragm wide open.


*4 Limited up to FN 22, no compliance with FN 26.5. *5 Analyzer and polarizer are recommended for usage with MPLFLN1.25X and 2.5X.

*5 Analyzer and polarizer are recommended for usage with MPLFLN1.25X and 2.5 *6 BD: Brightfield/darkfield objectives.

*7 Slight vignetting may occur in the periphery of the field when MPLN-BD series objectives are used with high-intensity light sources, such as mercury and xenon, for darkfield observation.

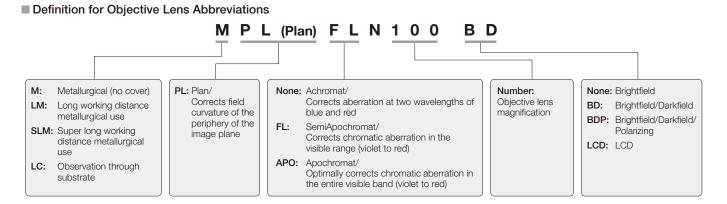

GX53 System Diagram

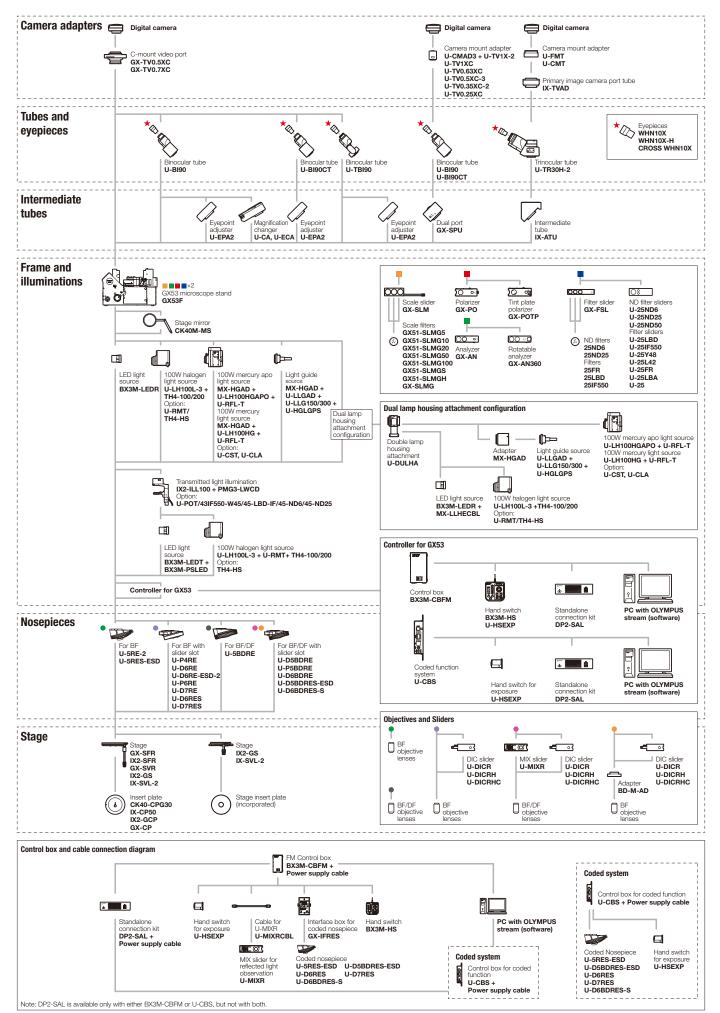


Specifications

| | | GX53 |
|------------------|-------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Optical system | | UIS2 optical system (infinity-corrected) |
| | Reflected light illumination | Manual brightfield/darkfield selection by mirror unit Manual field stop/aperture stop switch with centering Light source: White LED (with Light Intensity Manager) /12 V, 100 W halogen lamp/100 W mercury lamp/light guide source Observation mode: brightfield, darkfield, differential interface contrast (DIC)*1, simple polarizing*1, MIX observation (4 directional darkfield)*2 *1 Slider for exclusive use of this observation is required. *2 MIX observation configuration is required. |
| | Transmitted light illumination (optional) | Stand for transmitted light (IX2-ILL100: with field stop) is required PMG3-LWCD: Condenser (NA 0.6, WD 12 mm) with aperture stop Light source: White LED (with Light Intensity Manager) 12 V, 100 W halogen lamp Observation modes: brightfield, simple polarizing |
| | Imprinting of scale | All ports reversed positions (up/down) from observation positions seen through the eyepiece |
| | Output front port (optional) | Camera and DP system (reversed image, special camera adapter for GX) |
| | Output side port (optional) | Camera, DP system (upright image) |
| Microscope frame | Electrical system | Reflected light illumination Built-in LED power supply for reflected light illumination Continuously-variable light intensity dial Input rating 5 V DC, 2.5 A (AC adapter 100–240 V, AC 0.4 A, 50 Hz/60 Hz) Transmitted light illumination (requires the optional BX3M-PSLED power supply) Continuously-variable light intensity dial by voltage Input rating 5 V DC, 2.5 A (AC adapter 100–240 V, AC 0.4 A, 50 Hz/60 Hz) External interface (requires the optional BX3M-CBFM control box) Coded nosepiece connector × 1 MIX Slider (U-MIXR) connector × 1 Handset (U-HSEXP) connector × 1 RS-232C connector × 1, USB 2.0 connector × 1 |
| | Focus | Rack and pinion with roller guide Manual, coarse and fine coaxial handle; focus stroke 9 mm (2 mm above and 7 mm below the stage surface) Fine handle stroke per rotation: 100 µm (min. scale: 1 µm) Coarse handle stroke per rotation: 7 mm With torque adjustment ring for coarse focusing With upper limit stopper for coarse focusing |
| Tubes | Widefield (FN 22) | Inverted: binocular (U-BI90, U-BI90CT), trinocular (U-TR30H-2), tilting binocular (U-TBI90) |
| Nosepiece | | Brightfield Holes: 4 to 7 pcs, Type: Manual/Coded, Centering: Enabled/Disabled Brighfield/darkfield Hole: 5 to 6 pcs, Type: Manual/Coded, Centering: Enabled/Disabled |
| Stage | | Right handle stage for GX (X/Y stroke: 50 × 50 mm, max. load 5 kg) Flexible right handle stage, left short handle stage (each X/Y stroke: 50 × 50 mm, max. load 1 kg) Gliding stage (max. load 1 kg) A set of teardrop and long hole types |
| Weight | | Approx. 25 kg (microscope frame 20 kg) |
| Environment | | Indoor use Ambient temperature: 5 to 40 °C (45 to 100 °F) Maximum relative humidity: 80% for temperatures up to 31 °C (88 °F) (without condensation) In case of over 31 °C (88 °F), the relative humidity is decreased linearly through 70% at 34 °C (93 °F), 60% at 37 °C (99 °F), and to 50% at 40 °C (104 °F). Pollution degree: 2 (in accordance with IEC60664-1) Installation/Overvoltage category: II (in accordance with IEC60664-1) Supply voltage fluctuation: ±10 % |

Dimensions

GX53

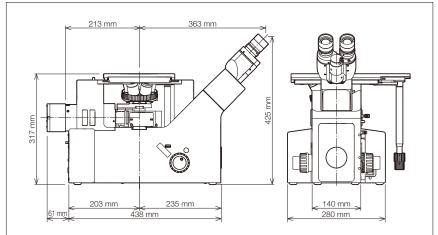

Olympus offers an extensive product line for materials science and industrial microscopy. Learn more about the LEXT 3D measuring laser microscope and DSX series digital microscopes on our website, www.olympus-ims.com.



NEW

LEXT 3D Measuring Laser Microscope

Power and speed are the hallmarks of the LEXT OLS5000 microscope. Owing to its high-resolution imaging and fast acquisition, the microscope delivers precise non-contact 3D observations and measurement of surface topography with impressive efficiency.



DSX Digital Microscope

The DSX series' advanced digital technology delivers surpassing image quality with operating simplicity, making it suitable for users of any experience level. The DSX microscope's intelligent interface is as easy-to-use as using a smartphone or tablet and backed by guaranteed accuracy and repeatability for 2D and 3D measurements.

· OLYMPUS CORPORATION is ISO14001 certified.

- · OLYMPUS CORPORATION is ISO9001 certified.

- This product is designed for use in industrial environments for the EMC performance. Using it in a residential environment may affect other equipment in the environment.
 All company and product names are registered trademarks and/or trademarks of their respective owners.
 Images on the PC monitors are simulated.
 Specifications and appearances are subject to change without any notice or obligation on the part of the manufactur
 Illumination devices for microscope have suggested lifetimes. Periodic inspections are required.
 Please visit our web site for details.

www.olympus-ims.com

